

D12 Lab Flotation Machine KBDA-24D Variable Frequency Drive Setup

WARNING! HIGH VOLTAGE -

Disconnect main power before making connections.

Terminals “L1” and “L2” are for single phase AC line input only. Jumper “J1” should be set to the 115V position for use with a 115V power supply. For use with a 230V power supply, move jumper ”J1” to the 230V position.

RECONDITIONING THE BUS CAPACITORS

If this drive has been in storage for over one year it is necessary to recondition the power supply bus capacitors. To recondition the bus capacitors, apply the AC Line, with the drive in the Stop Mode, for a minimum of one hour. Not following this procedure will cause the bus capacitors to fail.

(A40194) – Rev. A00 – 7/31/2013

